

OAKTON® Waterproof pHTestr 3 & 3+ Double Junction Instructions

Before you Begin:

Remove electrode cap. To condition electrode, immerse electrode in electrode storage solution, buffer or tap water for at least 30 minutes. DO NOT use de-ionized water.

Calibration:

Calibration should be done regularly, typically every day that the Testr is used. You can calibrate at up to three points (pH 4.01, 7.00, and 10.00).

1. Press ON/OFF button to switch unit on.
2. Dip electrode 1/2" to 1" into buffer (pH 4.01, 7.00, or 10.00).
3. Press CAL button to enter Calibrate (CAL) mode. 'CAL' flashes on the display. Then, a pH value close to the pH buffer value will flash repeatedly.
4. After at least 3 minutes press HOLD/CON to confirm calibration. The display will show 'CO' and then switch to the buffer value reading.
5. Repeat with other buffers if necessary. Rinse electrode in tap water before dipping into next buffer.

Calibration Troubleshooting:

Failure to press HOLD/CON to confirm calibration (step 4 above). Pressing the CAL button will resume measuring mode but will not enter the calibration value.

Insufficient sampling time. If the Testr does not have a long enough exposure to the buffer, a stable calibration point will not be reached. Wait at least 3 minutes before pressing HOLD/CON.

Failure to re-hydrate the electrode. A dry electrode will give fluctuating readings while it rehydrates in a buffer, causing errors.

Buffer Set Selection:

To choose USA (4.01, 7.00, 10.00) or NIST (4.01, 6.86, 9.18) buffers:

1. With Testr off, hold HOLD/CON key while pressing ON/OFF key.
2. Press HOLD/CON key again.
3. Press CAL key to toggle between USA and NIST sets.
4. Press HOLD/CON to confirm.

Displaying Temperature (°C) for pHTestr 3+ only:

1. With Testr on, press HOLD/CON key. Then press CAL key. Testr will display °C.
2. Press HOLD/CON key again to return to pH mode.

pH Testing:

1. Remove cap from the electrode and press the ON/OFF button to switch Testr on.
2. Dip the electrode 1/2" to 1" into the test solution. Stir once and let the reading stabilize.
3. Note the pH or press HOLD/ CON button to freeze the reading. Press HOLD/CON again to release the reading.
4. Press ON/OFF to turn off Testr. If you do not press a button for 8.5 minutes the Testr will automatically shut off to conserve batteries.


Instrument Maintenance:

- Rinse the electrode with tap water or electrode storage solution after each measurement.
- Periodic soaks in warm pH 4.01 buffer will help remove contaminants.
- If possible, keep a small piece of paper or sponge in the electrode cap—moistened with clean water or electrode storage solution (NOT de-ionized water)—and close the cap over the electrode.

When you need a new electrode, see "Electrode Replacement" at right.

Changing Batteries:

1. Open battery compartment lid (with attached lanyard loop).
2. Remove old batteries; replace with fresh ones. Note polarity (shown in battery compartment and in picture at right).

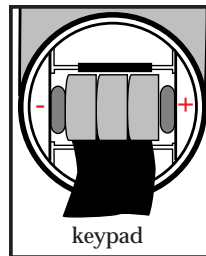

Error Messages:

bAt Weak batteries—replace

Er1 Wrong or bad buffer value, or electrode is failing.

or Ur Over range/under range signal, or electrode is not in contact with solution, or electrode is failing.

Specifications

	WP pHTestr 3	WP pHTestr 3+
pH		
Range	0.00 to 14.99 pH	
Resolution	0.01 pH	
Accuracy	±0.01 pH	
Calibration	up to 3 points (4.01; 7.00; 10.00 or 4.01; 6.86; 9.18)	
Temperature		
Range	—	0 to 50°C
Resolution	—	1°C
Accuracy	—	±1°C

ATC: 32 to 122°F/0 to 50°C

Operating temperature: 32 to 122°F/0 to 50°C

Special Functions: On/Off; HOLD; CA (Calibrate); CO (Confirm display); auto-buffer recognition; auto-shutoff after 8.5 min. nonuse

Dimensions: 6.5" L x 1.5" Dia. (165 x 38 mm)


Weight: 3.25 oz (90 gms)

Power: Three 1.5 V batteries (included). 250 hours continuous use.


Electrode replacement:

You can replace the electrode module at the fraction of the cost of a new Testr. When the Testr fails to calibrate, gives fluctuating readings in buffers, shows error messages 'E2' or 'OR' in a buffer, and the procedures in the Maintenance section do not help, you need to change the electrode.

1. With dry hands, grip the ribbed Testr collar with electrode facing you. Twist the collar counter clockwise. (see diagram A). Save the ribbed Testr collar and O-ring for later use.

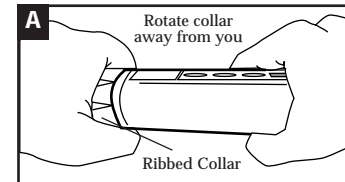

2. Pull the old electrode module away from the Testr.
3. Align the four tabs on the new module so they match the four slots on the testr. (see diagram B).

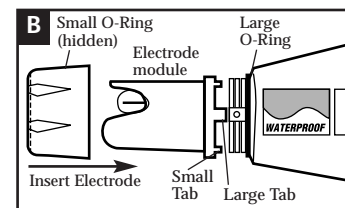

4. Gently push the module onto the slots to seat it in position. Push the smaller O-ring fully onto the new electrode module. Push the collar over the module and thread it into place by firmly twisting clockwise.

Warranty:

The waterproof pHTestr 3 meter body is warranted against defects in materials and workmanship for a period of 12 months from the date of purchase; the electrode module is warranted for a period of 6 months from the date of purchase. If repair, adjustment or replacement is necessary and has not been the result of abuse or misuse within the 6 month period, please return the Testr—freight pre-paid—and correction will be made without charge. Out of warranty products will be repaired on a charge basis.

Return of Items:

Authorization must be obtained from your OAKTON Distributor before returning items for any reason. When applying for authorization, please include information regarding the reason the item(s) are to be returned.

Note: We reserve the right to make improvements in design, construction and appearance of products without notice. Prices are subject to change without notice.